


# GlutenTox•Pro Surface Fast


## Quick Guide

**REF** KIT3003 (KT-6406)

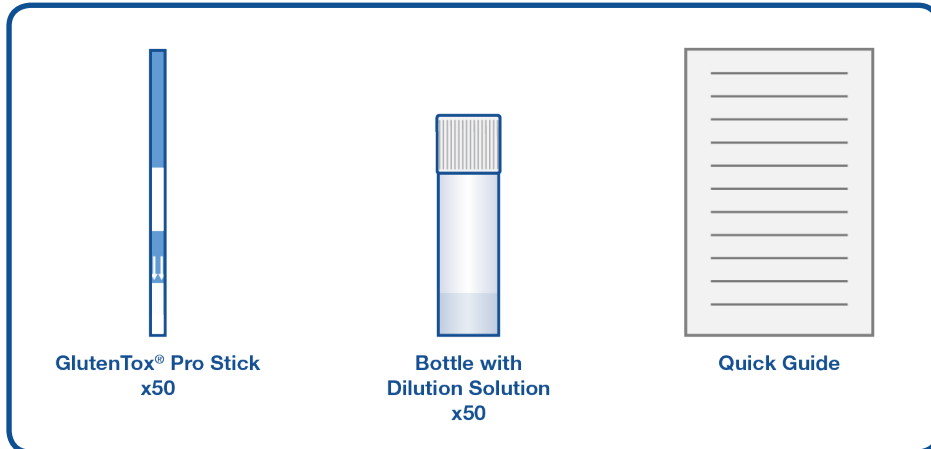

### Test procedure

1. Rub the white, cotton wool side of the stick against five areas of 1 cm<sup>2</sup>/ 0.16 in<sup>2</sup> or along a rectangle of 5 x 1 cm/0.16 x 1.97 in. Choose 5 representative areas distributed along the surface.
2. Open the dilution solution bottle and put the white end of the stick in contact with the liquid.
3. Let it stand in the bottle and wait 10 minutes. If a red line appears at 4 minutes, the working surface has a content in gluten equivalent to > 10 ppm.

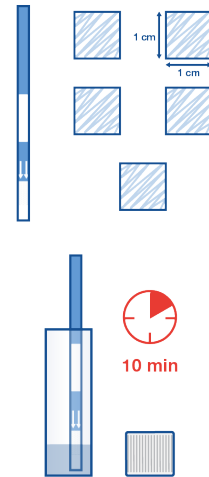

### Results interpretation

1. Final reading at 10 minutes

2. Initial reading at 4 minutes



**Negative**  
Clean



**Positive**  
Contaminated



**Invalid**



Control line  
Test line

**Positive**  
>10 ppm gluten