



Anton Paar

DMA 35

Standard Operating Procedure



Installation

Installation | Instalación | Installation | Instalação |
Kurulum | インストール | 安装



Water Check

Wasser-Check | Análisis del agua | Vérification à l'eau |
Verificação com água | Su Doğrulaması | 水によるチェック |
水検



Setting the Measuring Method

Auswahl der Messmethode | Configuración del método de
medición | Réglage de la méthode de mesure | Definir a
método de medição | Ölçüm Metodunu Ayarlama |
測定メソッドの設定 | 设置测量方法



Sample Identification

Probenidentifikation | Identificación de la muestra |
identification des échantillons | Identificação da amostra |
Örnek Tanımlama | サンプルの識別 | 样品标识



Degassing

Entgasen | Desgasificación | Dégazage | Desgaseificação |
Degaze etmek | 脱ガス | 脱気



6

Filling

Füllen | Llenado | Remplissage | Enchimento |
Doldurma | 充填 | 进样



7

Measurement

Messung | Mediciones | Measure | Medição |
Ölçüm | 測定 | 測量



8

Viewing Measured Data

Ansehen der Messdaten | Observación de los datos
medidos | Visualisation des données mesurées |
Visualizando os dados medidos | Ölçülen veriyi
görüntüleme | 測定結果の閲覧 | 查看测量数据



9

Handling Measured Data

Handhabung der Messdaten | El manejo de datos
de medición | Manipulation des mesurées données |
Manipulação de dados medidos | Ölçüm Sonuçlarını
Değerlendirme | 測定データの処理 | 处理测量数据








10


Cleaning

Reinigung | Limpieza | Nettoyage | Limpeza |
Temizleme | 洗净 | 清洗


Installation

1 ● ● ● ● ● ● ● ● ● ●

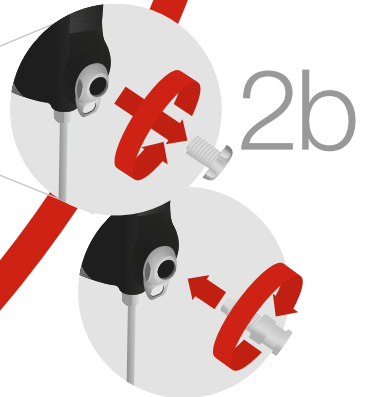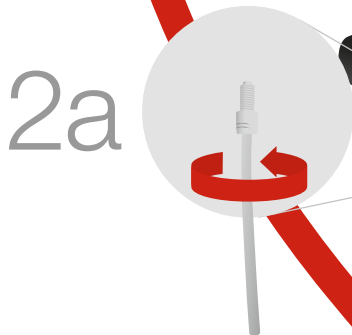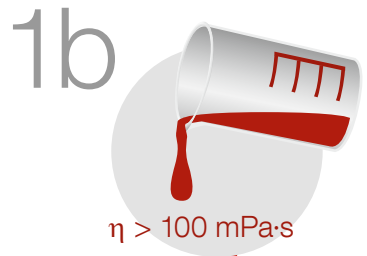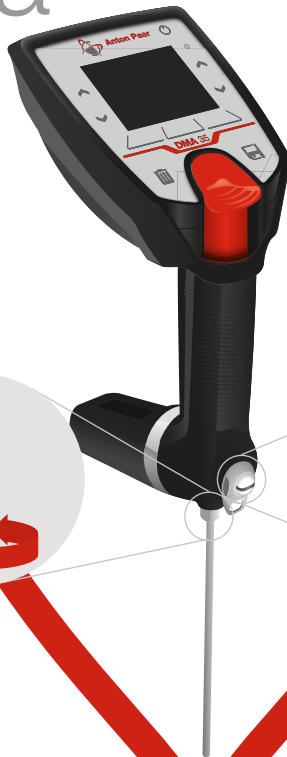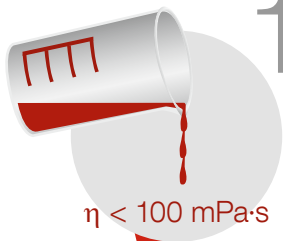


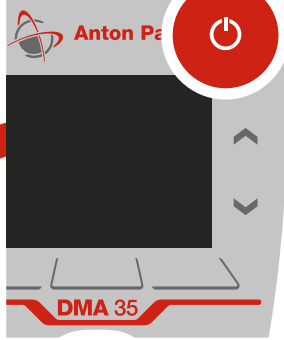

4

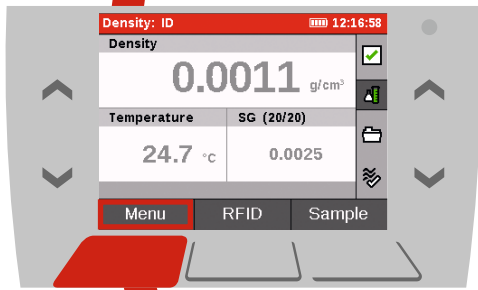


5

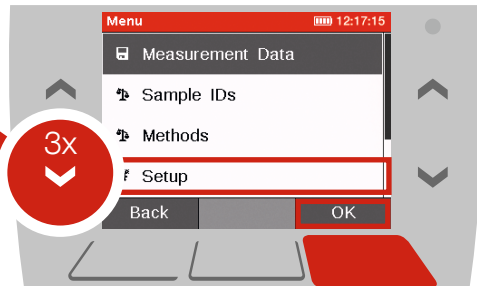

6


Water Check

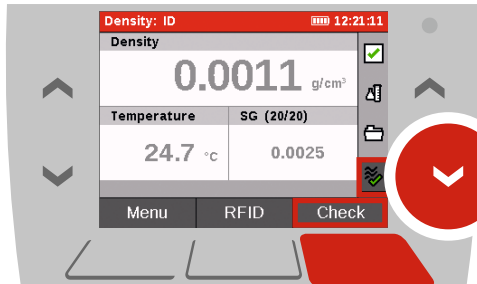

3



4

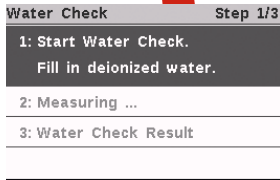

1

2



5






6





Deionized water*

* Entionisiertes Wasser | Agua desionizada | Eau déionisée | Água deionizada | İyonuz su | Uygun temizleme sıvısı | 脱イオン水 | 去离子水


7

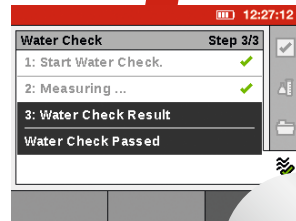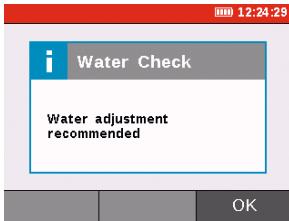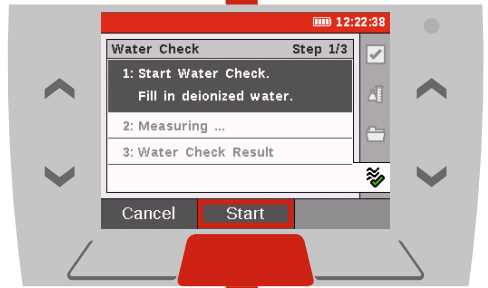

Setting the Measuring Method

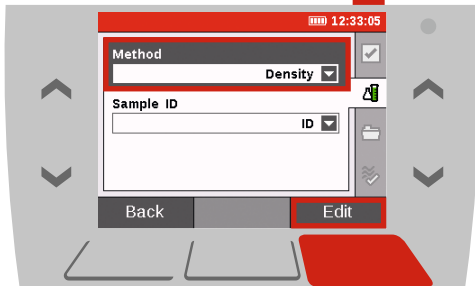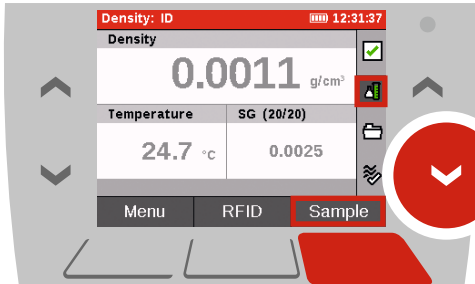


Sample Identification

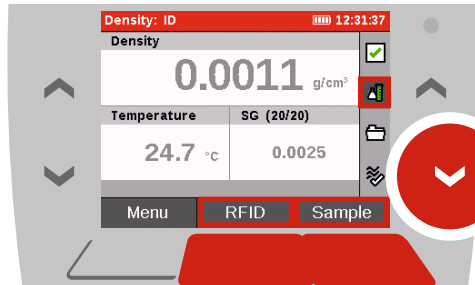

1a

1b


2a

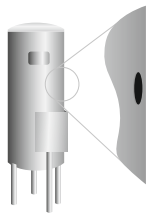

2b



RFID-Setup*:
Manual

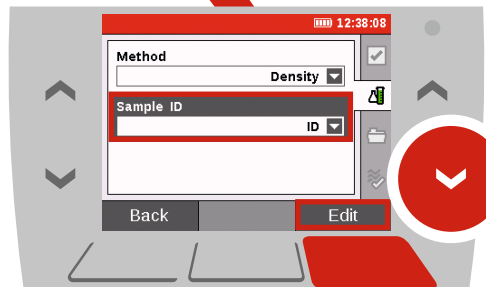


3a

4a


3b

Reading Tag...

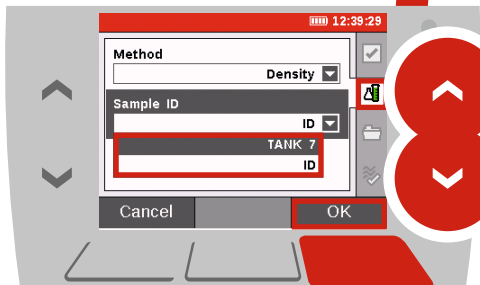

5a


6a

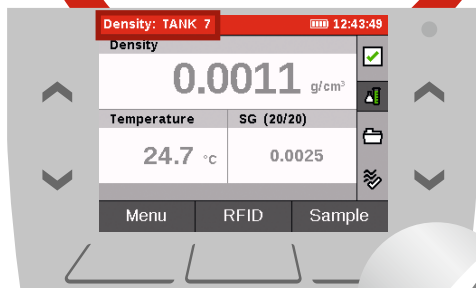

* Conexión RFID: Manual | Paramétrage RFID: Manuel Configuração do RFID: Manual | RFID-Ayarı: Kullanım Kılavuzu | RFID設定:マニュアル | RFID設置: 手冊

Degassing

5 ●●●●●●●●●●

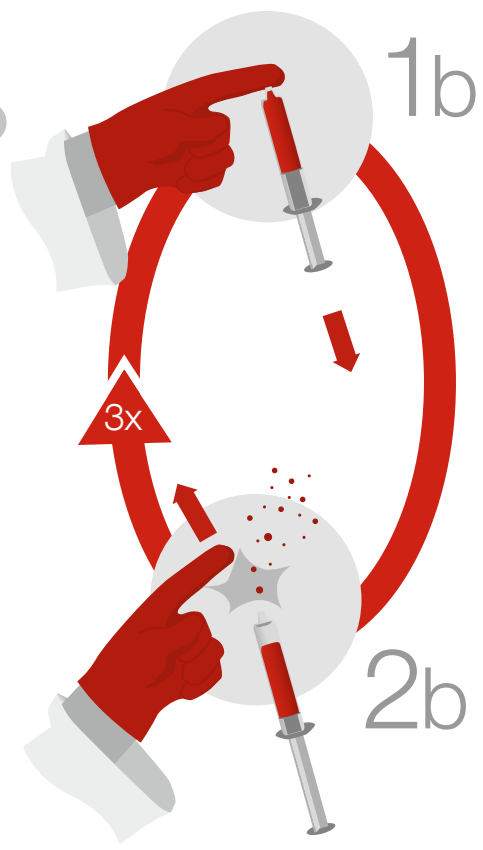

Filling




1a



2a




3a

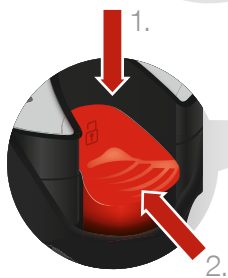

1b

Measurement

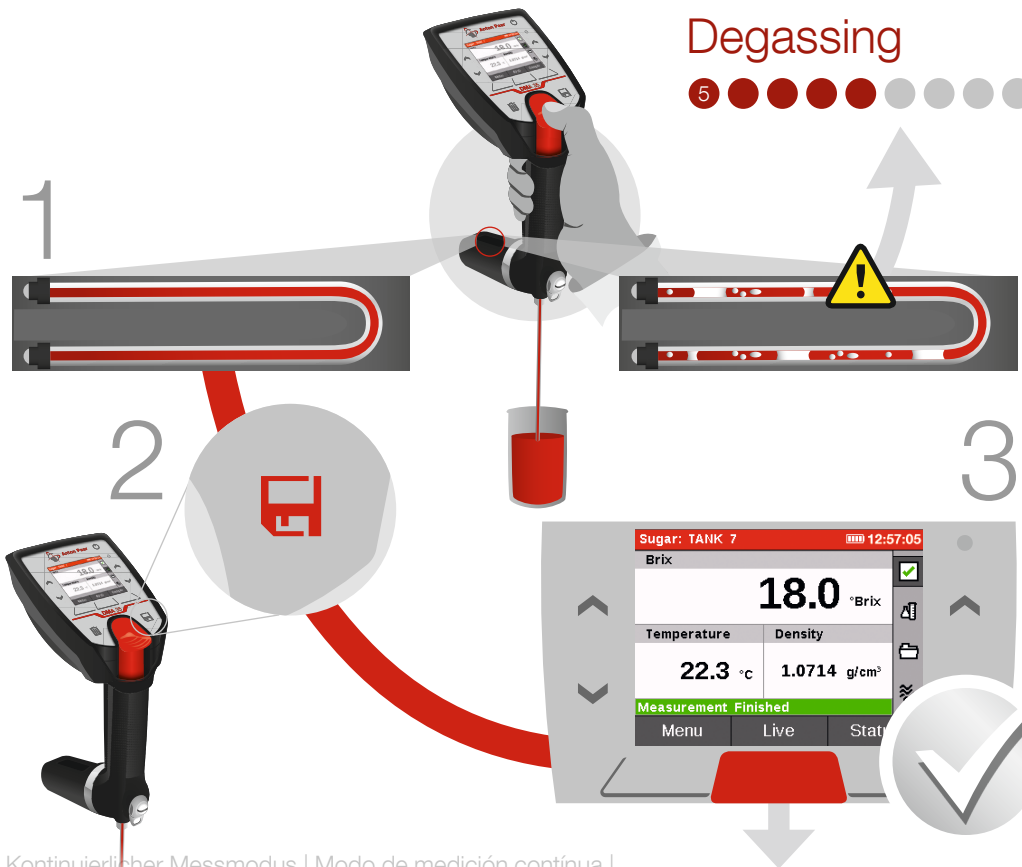


* Kontinuierlicher Messmodus | Modo de medición continuo |
Mode de mesure continu | Modo de medição contínuo |
Aralıksız Ölçüm Modu | 測定モードへ | 连续测量模式


Continuous
Measuring Mode*

Viewing Measured Data

8

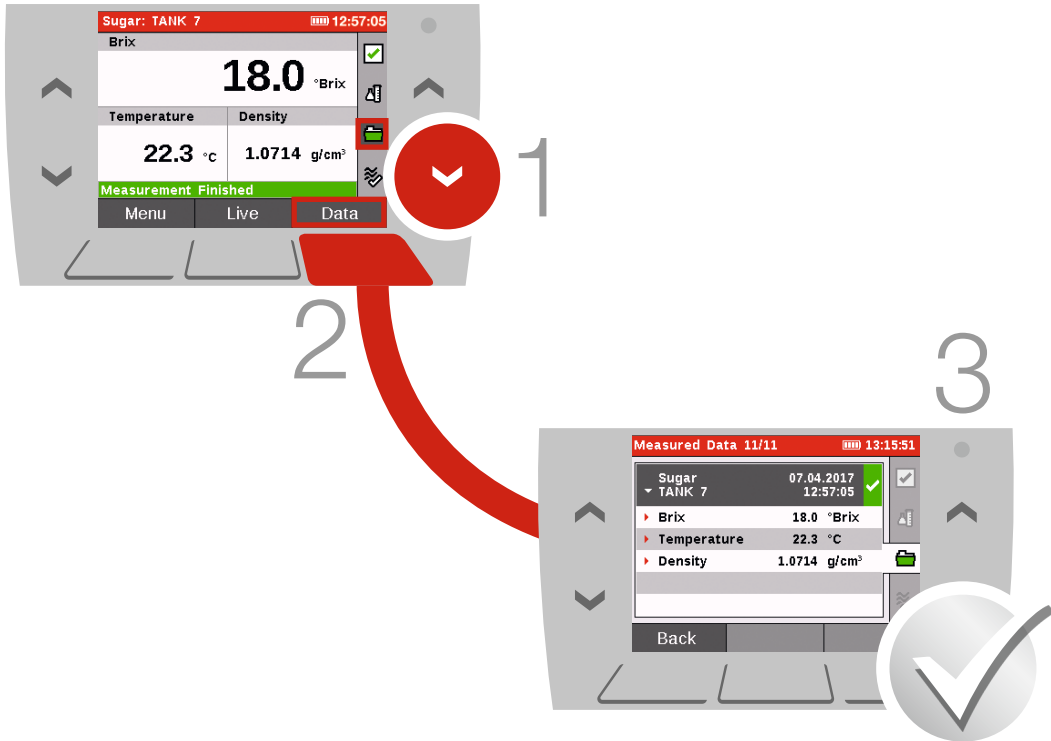

Handling Measured Data

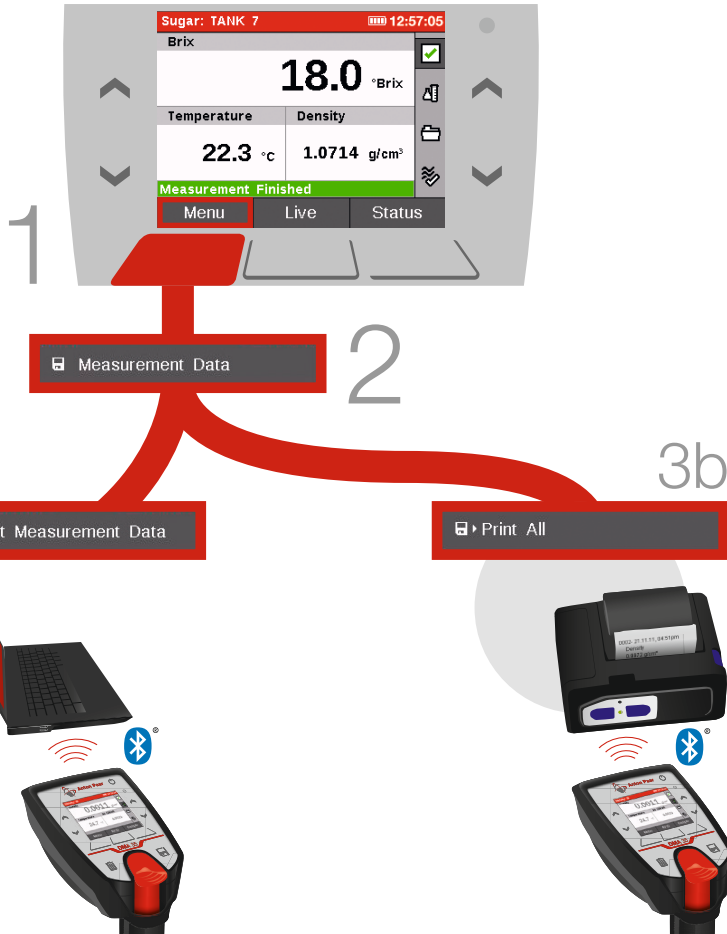

Cleaning

10 ● ● ● ● ● ● ● ● ● ●



* Geeignetes Lösungsmittel | Solvente adecuado | solvant approprié | Solvente adequado | Uygun temizleme sıvısı | 適切な溶媒 | 合适的清洗溶剂

Notes

© 2019 Anton Paar GmbH | All rights reserved.
Specifications subject to change without notice.
E281B003ML-D

www.anton-paar.com