



ONE SYSTEM BASIC


SINGLE-USE WIPE DISPENSER
FOR USE WITH **Rely+On®**
Virkon® DISINFECTANT

PRODUCT INFORMATION:

- Supplied in **packs of 4** dispensers
- **120 wipes** per dispenser
- **2.5l** fill capacity
- Dispenser dimensions: **17.5cm x 36cm**
- For use with **Rely+On® Virkon®** disinfectant
- Active for use in **15 minutes**
- Up to **5 days*** of standing time
- Filled with **DESOTEX** non-woven wipes
- **Compact**, flexible packaging
- **High stability** (Free-standing)
- **Easy handling** and fill-level control
- Tear-resistant, **lint-free**, non-woven

MIXING AND USAGE DIRECTIONS

- Add 25g **Rely+On® Virkon®** Powder (or 5 x 5g if using tablets) to 2.5l of water.
- Stir until **FULLY** dissolved.
- Add to **ONE SYSTEM BASIC** container as directed opposite
- Wipe surfaces to be disinfected, **leave for 10 minutes** (contact time), then wipe dry or rinse** with clean water.

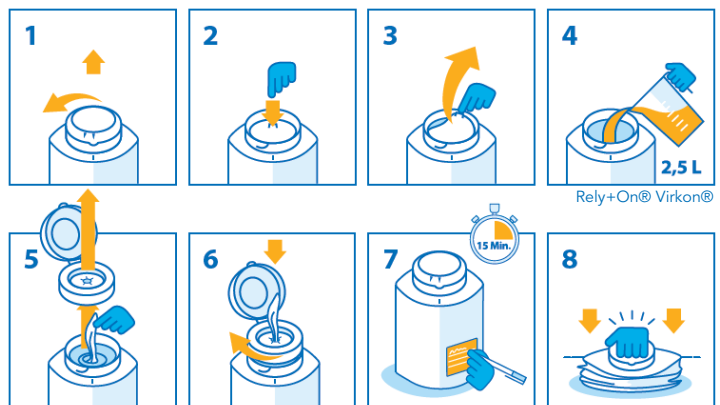

- Always store out of direct sunlight
- Rely+On® Virkon® 1% solutions are effective for up to 5 days*

In all instances, perform a risk assessment, then wear the appropriate level of personal protective equipment (when preparing and using disinfectant solutions).

* Valid when used with the supplied brand of DESOTEX wipes

** Rinse any sensitive surfaces with clean water after the disinfectant application contact time (refer to Rely+On® Virkon® materials compatibility guide).

Not for use on any medical devices. Contact Day-Impex for advice on use with suitable DESOTEX dry wipe systems.

Use biocides safely. Always read the label and product information before use.